EN - English	1
FR - Français	7
ES - Español	14
IT - Italiano	21
DE - Deutsch	28
PT - Português	35
NL - Nederlands	42
NO - Norsk	46
DA - Dansk	50
SV - Svenska	54
FI - Suomi	58
HU - Magyar	62
PL - Polski	66

European Catalogue Numbers:

Euro Plug: 444-2900

UK Plug: 444-2901

Swiss Plug: 444-2902

Table of ConTenTs

Package contents	1
Warranty	1
Installation.....	2
Maintenance & servicing.....	2
Intended use	2
Equipment disposal	2
Environmental conditions.....	2
Safety instructions.....	3
Warning	3
Standard & regulations	3
Specifications.....	4
Operating instructions	5
Technical service.....	6
Troubleshooting	6
EC Declaration of Conformity.....	70-71

PaCkage ConTenTs

Orbital Shaker
 Non-skid rubber mat,
 234cm detachable power cord
 Instruction manual

WarranTy

VWR International warrants that this product will be free from defects in material and workmanship for a period of two (2) years from date of purchase. If a defect is present, VWR will, at its option, repair, replace, or refund the purchase price of this product at no charge to you, provided it is returned during the warranty period. This warranty does not apply if the product has been damaged by accident, abuse, misuse, or misapplication, or from ordinary wear and tear. For your protection, items being returned must be insured against possible damage or loss. This warranty shall be limited to the replacement of defective products. IT IS EXPRESSLY AGREED THAT THIS WARRANTY WILL BE IN LIEU OF ALL WARRANTIES OF FITNESS AND IN LIEU OF THE WARRANTY OF MERCHANTABILITY.

Installation

Upon receiving the VWR Orbital, check to ensure that no damage has occurred during shipment. It is important that any damage that occurred in transport is detected at the time of unpacking. If you do find such damage the carrier must be notified immediately. After unpacking, place the shaker on a level bench or table, away from explosive vapors. Ensure that the surface on which the unit is placed will withstand typical heat produced by the unit and place the unit a minimum of six (6) inches from vertical surfaces. Always place the unit on a sturdy work surface. The Orbital Shaker is supplied with a power cord that is inserted into the IEC connector on the back of the unit first, then it can be plugged into a properly grounded outlet. The 230V unit plugs into a 230 volt, 50/60 Hz source.

Maintenance & Servicing

The Orbital Shaker is built for long, trouble-free, dependable service. No lubrication or other technical user maintenance is required. It needs no user maintenance beyond keeping the surfaces clean. However at least every three (3) months you should:

- Unplug the unit.
- Remove any accumulated dirt from the base and tray.
- Check all accessible items to make sure they are properly tightened.

The unit should be given the care normally required for any electrical appliance. Avoid wetting or unnecessary exposure to fumes. **DO NOT** use a cleaning agent or solvent on the front panel which is abrasive or harmful to plastics, nor one which is flammable. Always ensure the power is disconnected from the unit prior to any cleaning. If the unit ever requires service, contact your VWR representative.

Intended Use

These Shakers are intended for general laboratory use.

Equipment

Disposal

This equipment must not be disposed of with unsorted waste. It is your responsibility to correctly dispose of the equipment at life-cycle-end by handing it over to an authorized facility for separate collection and recycling. It is also your responsibility to decontaminate the equipment in case of biological, chemical and/or radiological contamination, so as to protect the persons involved in the disposal and recycling of the equipment from health hazards.

For more information about where you can drop off your waste of equipment, please contact your local dealer from whom you originally purchased this equipment. By doing so, you will help to conserve natural and environmental resources and you will ensure that your equipment is recycled in a manner that protects human health.

Environmental Conditions - Standard

Operating Conditions: Indoor use only.

* For use in CO₂ environments, incubators, or cold rooms.

Temperature:	0 to 40°C
Humidity:	maximum 80% relative humidity, non-condensing
Altitude:	0 to 2000 M above sea level

Non-Operating Storage:

Temperature:	-20 to 65°C
Humidity:	maximum 80% relative humidity, non-condensing

Installation Category II and Pollution Degree 2 in accordance with IEC 664.

SafeTy InsTrUCTIONS

Please read the entire instruction manual before operating the Orbital Shaker.

WARNING! DO NOT use the Orbital Shaker in a hazardous atmosphere or with hazardous materials for which the unit was not designed. Also, the user should be aware that the protection provided by the equipment may be impaired if used with accessories not provided or recommended by the manufacturer, or used in a manner not specified by the manufacturer.

Always operate unit on a level surface for best performance and maximum safety.

DO NOT lift the unit by the tray.

CAUTION! To avoid electrical shock, completely cut off power to the unit by disconnecting the power cord from the unit or unplug from the wall outlet. Disconnect unit from the power supply prior to maintenance and servicing.

Spills should be removed promptly. **DO NOT** immerse the unit for cleaning.

DO NOT operate the unit if it shows signs of electrical or mechanical damage.

Earth Ground - Protective Conductor Terminal

Alternating Current

Pinch Point - Keep fingers clear during operation.

Standards

RegUlaTions

VWR international hereby declares under its sole responsibility that the products comply with the requirements of the following directives and associated standards:

Associated EU directives:

EMC directive	2014/30/EU
LVD directive	2014/35/EU
ROHS directive	2011/65/EU
WEEE directive	2012/19/EU

Safety standards:

IEC 61010-1
IEC 61010-2-051
EN 61010-1
EN 61010-2-051
EN 50581

EMC standards:

EN 61326-1

Overall dimensions LxWxH	27,9 x 43,2 x 10,2cm
Tray dimensions LxW	22,2 x 29,8cm
Electrical	230 volts, 2,5 amps, 25 watts
Fuses	5mm x 20mm, 5 amp quick acting
Speed Range	40 to 300 rpm
Orbit	15mm
Capacity	3,6kg
Ship Weight	11,3kg
European Catalogue Numbers:	
	Euro plug: 444-2900
	UK plug: 444-2901
	Swiss plug: 444-2902

OPERATING INSTRUCTIONS - Standard Orbital Shaker

The Standard Orbital Shaker is used for general laboratory shaking needs.

1. Getting ready:

a. The speed knob should be at the extreme counter-clockwise position or at the #1 on the dial.

b. Make sure the rocker switch is in the standby position.

c. Plug the cord into a properly grounded outlet.

2. Setting speed:

a. To run push rocker to the on position. Set speed knob to desired setting and if necessary. Unit will run until you move the rocker switch to the off position. The microprocessor speed control slowly ramps to set speed to avoid splashing.

3. Turning unit off:

a. To stop shaking function, turn the speed knob to the extreme counter-clockwise position and push the rocker switch to the standby position. The shaker should be kept in the standby position when not in use. To completely cut off power to the unit, disconnect the power cord from the unit or unplug from the wall outlet.

OPERATING TIPS

Centering your sample and even weight distribution on the tray helps with balance and stability.

The shaker will automatically restart after a power interruption.

TeChnIcal SerVICe

For more information or technical assistance contact your local VWR representative or visit www.vwr.com.

Web resources: Visit VWR's website at www.vwr.com for:

- Complete technical service contact information
- Access to VWR's Online Catalogue, and information about accessories and related products
- Additional product information and special offers

TroUbleShooTIng

Problem	Cause	Solution
Unit fails to power on	Missing or blown fuse	Add or replace fuse as necessary. If problem persists, please contact your VWR representative for repair.
Unit does not shake	Mechanical obstruction Drive system failure Ceased bearing Drive belt broken	Remove mechanical obstruction. If problem persists, the reason may be the drive system. Please contact your VWR representative for repair.
Unit is excessively noisy	Rattling or ticking sounds may indicate a loose screw on the tray	Ensure that tray is secured tightly. If problem persists, please contact your VWR representative for repair.
Unit stops running after initial start-up	Maximum load exceeded	Ensure the load is evenly distributed and does not exceed the maximum load capacity for the unit. If problem persists, please contact your VWR representative for repair.

EN - English	1
FR - Français	7
ES - Español	14
IT - Italiano	21
DE - Deutsch	28
PT - Português	35
NL - Nederlands	42
NO - Norsk	46
DA - Dansk	50
SV - Svenska	54
FI - Suomi	58
HU - Magyar	62
PL - Polski	66

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The undersigned states that this declaration is issued under the sole responsibility of the manufacturer, VWR International bvba, and that the object(s) of the declaration described in annex 1 are in conformity with the relevant Union harmonization legislation.

In the event of unauthorized modification of any of the products listed in the annex 1, this declaration becomes invalid.

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Document Number: 14871.2017.07

Place & date of issue: Leuven, 3/03/2017

Signed for and on behalf of VWR International bvba:

Yves Van Damme
European Regulatory Affairs Manager

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India

VWR Lab Products Private Limited
No.139, BDA Industrial Suburb,
6th Main, Tumkur Road, Peenya Post,
Bangalore, India – 560058
Tel.: +91-80-28078400
Email: vwr_india@vwr.com

Ireland / Northern Ireland

VWR International Ltd / VWR International
(Northern Ireland) Ltd
Orion Business Campus
Northwest Business Park
Ballycoolin
Dublin 15
Tel.: +353 01 88 22 222
Email sales.ie@vwr.com

Italy

VWR International S.r.l.
Via San Giusto 85
20153 Milano (MI)
Tel.: +39 02-3320311
Email: info.it@vwr.com

The Netherlands

VWR International B.V.
Postbus 8198
1005 AD Amsterdam
Tel.: +31 020 4808 400
Email: info.nl@vwr.com

Norway

VWR International AS
Haavard Martinsens vei 30
0978 Oslo
Tel.: +47 22 90 00 00
Email: info.no@vwr.com

Poland

VWR International Sp. z o.o.
Limbowa 5
80-175 Gdansk
Tel.: +48 058 32 38 200
Email: info.pl@vwr.com

Portugal

VWR International –
Material de Laboratório, Lda
Centro Empresarial de Alfragide
Rua da Indústria, nº 6
2610-088 Alfragide
Tel.: +351 21 3600 770
Email: info.pt@vwr.com

Singapore

VWR Singapore Pte Ltd
18 Gul Drive
Singapore 629468
Tel: +65 6505 0760
Email: sales.sg@vwr.com

Spain

VWR International Eurolab S.L.
C/ Tecnologia 5-17
A-7 Llinars Park
08450 - Llinars del Vallès
Barcelona
Tel.: +34 902 222 897
Email: info.es@vwr.com

Sweden

VWR International AB
Fagerstagatan 18a
163 94 Stockholm
Tel.: +46 08 621 34 00
Email: info.se@vwr.com

Switzerland

VWR International AG
Lerzenstrasse 16/18
8953 Dietikon
Tel.: +41 044 745 13 13
Email: info.ch@vwr.com

UK

VWR International Ltd
Customer Service Centre
Hunter Boulevard
Magna Park
Lutterworth
Leicestershire
LE17 4XN
Tel.: +44 (0) 800 22 33 44
Email: uksales@vwr.com

Legal Manufacturer:

VWR International bvba • Researchpark Haasrode 2020 • Geldenaaksebaan 464 • B-3001 Leuven • + 32 16 385011 • <http://be.vwr.com>
715091-00 (Rev 9)