# accu-jet®-Pipettierhelfer accu-jet® Pipette Controller

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Read carefully before use!

## Safety Instructions

This instrument may be involved with hazardous materials. This Operating Manual does not purport to address all of the safety problems associated with its use. It is the responsibility of whoever uses this instrument to consult and establish appropriate safety and health practices and to determine the applicability of regulatory limitations prior to use.

1. Follow general lab safety precautions and regulations, e.g., wear protective clothing, goggles and gloves, etc.
2. Observe all precautions from the Material Safety Data Sheet (MSDS) for any reagent you use with this product.
3. Every user must read and understand this Operating Manual before using the instrument.
4. Never use the instrument in an atmosphere that might be explosive.
5. Always work in a manner which neither endangers the user, nor any other person.
6. Use the instrument only for pipetting, and within the recommended limits of its chemical resistance and mechanical properties.
7. Never use force when using this instrument.
8. In case of instrument malfunction, immediately stop pipetting. Clean and repair the instrument according

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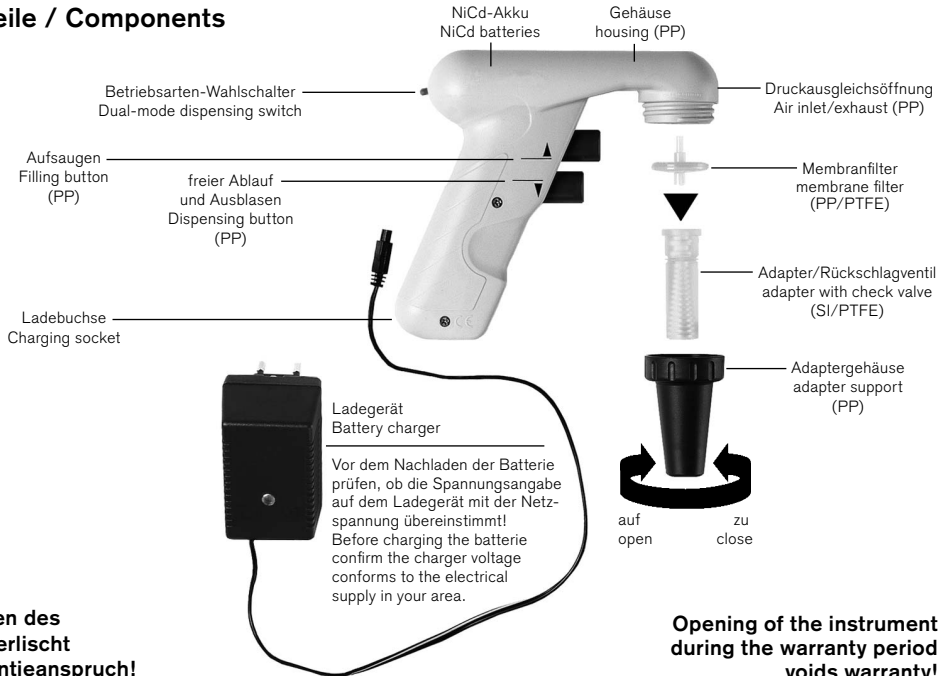
to the instructions before any further use of the instrument (➤ pages 13-18) or contact an authorized dealer or manufacturer.

9. Only use original manufacturer's accessories and spare parts.
10. Only the original recharger can be used for recharging of the NiCd batteries.
11. The recharger has to be protected against moisture and must only be used for this instrument.
12. Repairs other than battery changes are to be performed by trained and authorized service personnel only.
13. Opening of the instrument during the warranty period voids the warranty. If there is a failure during the warranty period, contact BrandTech for warranty service.
14. The instrument must only be used with original manufacturer's NiCd batteries.
15. Batteries must be completely discharged before disposal. Disposal should be performed in accordance with all governmental regulations.

### **Warning!**

Improper use of the instrument or the batteries such as short-circuiting, mechanical damage, overheating, may cause an explosion.

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## Storage conditions

Instrument and battery charger should be stored in a cool, dry place.

### **Storage temperature:**

- 20 °C to + 50 °C

## Starting operation

Before starting to operate, check to confirm the battery charger is of the appropriate voltage for your area and recharge the batteries, ➤ pages 16-17.

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## Instrument application

The instrument is designed to assist the filling and dispensing of graduated and volumetric pipettes of glass or plastic used for measuring liquids in the volume range of 0.1 mL to 100 mL.

It is not designed for Pasteur pipettes.

If the instrument is used correctly, the pipetted liquid will only contact the pipette.

## Restrictions of use

The instrument must not be used with liquids whose vapors are incompatible with silicone, PTFE, PP, EPDM and PF (phenol-formaldehyde resins).

### Warning!

This instrument must not be used or recharged in an explosive atmosphere. Highly flammable liquids (e.g., ether, acetone and other liquids with a flash point below 0 °C) must not be pipetted.

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## Restrictions of temperature

### Working- and recharging temperature:

+ 10 °C to + 35 °C

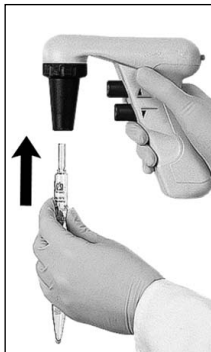
## Pipetting

### Attach the pipette

Prior to attaching the pipette, inspect the pipette's upper end for damage. A chipped pipette may damage the silicone pipette adapter. Hold the pipette as near to its upper end as possible, and carefully insert it into the adapter until it fits tightly.

#### **Attention!**

Be sure the pipette fits tightly into the adapter. Excessive force can cause the pipette to break. Use extra caution when inserting thin pipettes.

Once the pipette has been securely attached, always hold the instrument in a vertical position, tip down.

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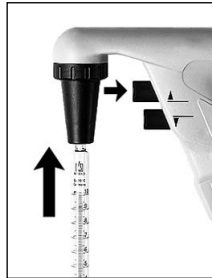
## Fill the pipette

Immerse the pipette tip into the liquid. Slowly press the upper pipetting button. The speed of aspiration is dependent on the amount the pipetting button is depressed. Fill the pipette so that the meniscus is slightly above the mark desired.

### Attention!

Take care to avoid aspiration of liquid into the instrument, which can impair instrument performance.  
If liquid is aspirated, clean the pipette adapter/check valve, and replace the membrane filter, if necessary (➤ pages 13-16).

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## Adjust the volume

Use suitable lint-free tissue to wipe the pipette tip. Slowly press the lower button until the meniscus is adjusted exactly to the desired volume.

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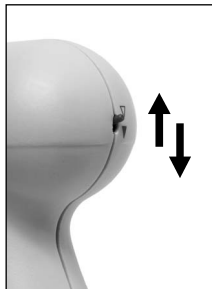
## Dispense the pipette

Select the appropriate dispensing mode

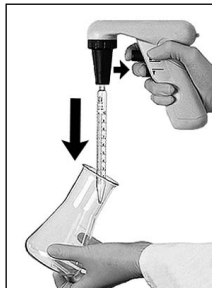
Switch up:

### ▽ Gravity-driven delivery

In this mode, the instrument allows the pipette to retain the volume of liquid prescribed for the accuracy of gravity driven-delivery pipettes.

Switch down:

### ▼ Blow-out (delivery with motor power)

## For both modes:

Slowly press the lower pipetting button. The speed of dispensing is dependent upon the amount the dispensing button is depressed.

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## Instrument storage

Attach the instrument to the wall support, as shown, for storage.

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### Note:

To mount the wall support, first clean the desired mounting area with a 50 % isopropyl alcohol / 50% deionized water mixture. (The surface of the area must be smooth). Remove the protective film from the snap tapes and tightly press the wall support to the surface. The snap tapes allow the wall support to be easily removed for cleaning. The wall support can also be mounted with two screws.

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## Cleaning

When necessary, wipe the polypropylene housing with a damp cloth.

When properly used, the instrument is maintenance-free.

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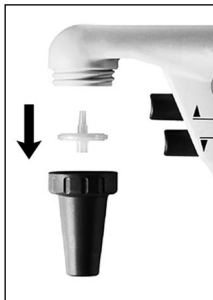
## Cleaning the check valve and replacing the filter

### Attention!

Always wear protective gloves and safety goggles!

In the event of pipette overfilling, or if the performance of the accu-jet® is impaired, observe the following instructions:

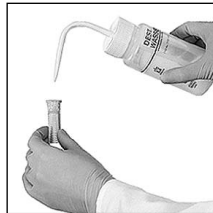
1. Remove pipette.
2. Gently unscrew the adapter support.
3. Remove the membrane filter.

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4. Remove the adapter from the adapter support by pushing it upwards. Do not disassemble the silicone pipette adapter!

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5. To clean the check valve, rinse the adapter through, using a wash bottle filled with water or other cleaning solution compatible with silicone and PTFE. Blow out the liquid, and allow drain completely.

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7. To insert the filter, the larger diameter connection must be facing the pipette adapter, and the smaller diameter connection must face the instrument body.

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8. The pipette adapter, adapter support and membrane filter may be steam sterilized at 121 °C for 20 minutes.

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**Note:**

Not all membrane filters are suitable for use with the instrument. For proper performance, only use the genuine manufacturer's recommended replacement filters.

Each instrument is supplied with a 0.2 µm membrane filter. The membrane filter can be steam sterilized up to 5 times at 121 °C.

It is the user's responsibility to ensure proper sterilization.

Use of the 0.2 µm (cell culture) filter allows for pipetting to Class B tolerances. For class A accuracy, use of 3 µm filter is necessary.

The larger pore size allows for proper gravity delivery times.

(Ordering information 3 µm filter, > page 27).

9. Reassemble the instrument by reversing the steps above.
10. To test the instrument for leaks, see "Leak test" (> page 16).

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## Leak Test

When the meniscus has been set, liquid should not drip out of the pipette until the valves are activated. Should the pipette drip, disassemble the pipette adapter and reassemble it carefully.

## Recharging the batteries

One full charge of the NiCd batteries allows 8 hours of non-stop pipetting. Decreasing motor power indicates the need for recharging.

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### Warning!

Use original battery charger only!  
Use of a different charger could damage the battery and would void the warranty. Confirm the voltage of the charger matches the voltage of your electrical outlet. The power supplied to the charger should differ by no more than  $\pm 10\%$ .

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## Recharging procedure

1. **Temperature range for recharging:** + 10 °C to + 35 °C
2. A light-emitting diode (LED) on the charger will glow while the accu-jet® is recharging.
3. For the first six hours, the instrument will charge in a high-current mode. Afterwards, the charger switches to a pulsed charge for long term battery life (LED will flash; 5 sec. on, 20 sec. off).
4. After 16 hours the batteries will be fully charged.

### Note:

- The instrument can be used during recharging.
- If the LED does not light up when the accu-jet® is connected to the charger, it will not charge. The accu-jet® and charger should both be tested and, if necessary, repaired or replaced.
- If the accu-jet® does not charge and the LED is lit, the battery should be replaced.

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## Battery replacement

### Note:

Use only the original manufacturer's replacement battery (ordering information > page 28).

### Warning!

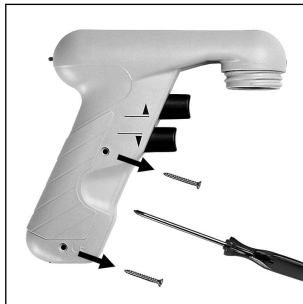
Use of the wrong batteries or improper use of the instrument (e.g., short-circuit, mechanical damage, overheating etc.) may cause the batteries to explode.

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

Lay the instrument on its left side  
(see photo). Use the point of a  
Phillips-head screwdriver to pierce  
the sealing wax covering the two  
screws holding the housing together.  
Remove the screws.

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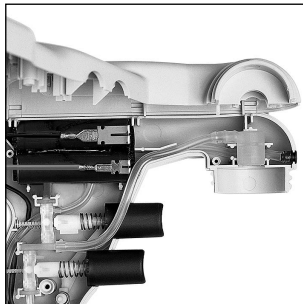Gently lift the right side housing and  
unfold the housing along the top axis  
(see photo).

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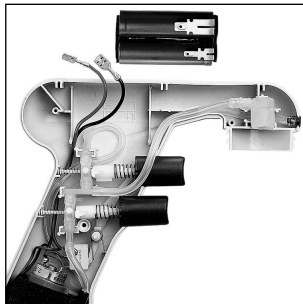
Gently slide the battery spade connectors apart and remove the battery.

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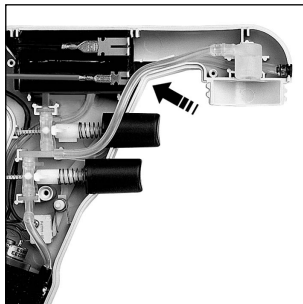
Insert the new battery and reconnect the spade connectors as far as they will go. Push the wires to the bottom of the housing.

**Ensure that the tubing is in the correct position so that it does not get pinched by the edge of the housing (see photo).**

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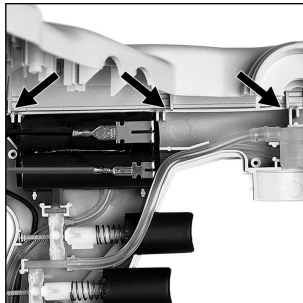Align the top of the housing. Make sure the three guide pins have the same placement as in the photo.

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When closing the housing it may be necessary to push the tubing inside the instrument. Use the end of your screwdriver to assist.

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Flip the dispensing mode switch to the lower position (powered dispensing). Test the pipetting buttons to check for motor function. Replace the screws and reattach the adapter support.

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## Disposal

NiCd-batteries must be completely discharged before disposal. Disposal must be carried out in accordance with governmental regulations. For disposal of the instrument, remove the batteries (➤ pages 18-20), and dispose of them separately.

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## Troubel Shooting

Trouble	Possible Cause	Action to be taken
Instrument does not aspirate properly	<ul style="list-style-type: none"> <li>- Check valve closed</li> <li>- Filter or check valve dirty</li> <li>- Battery discharged</li> </ul>	<ul style="list-style-type: none"> <li>- Push the switch to the lower position (blow out mode), and press the dispensing button (➤ p. 11)</li> <li>- Clean the check valve and replace the filter, if necessary (➤ p. 13-16)</li> <li>- Recharge the battery (➤ p. 16-17)</li> </ul>
Filling only possible after initial blow out, pressing the lower (dispensing) button	<ul style="list-style-type: none"> <li>- Filter or check valve dirty</li> </ul>	<ul style="list-style-type: none"> <li>- Clean check valve and replace the filter if necessary (➤ p. 13-16)</li> </ul>
Pipette drips	<ul style="list-style-type: none"> <li>- Filter installed upside down</li> <li>- Pipette or adapter damaged</li> </ul>	<ul style="list-style-type: none"> <li>- Reinstall filter (➤ p. 14)</li> <li>- Check pipette or adapter for damaged (➤ p. 13-16)</li> </ul>
Pipette not held properly	<ul style="list-style-type: none"> <li>- Adapter dirty or damaged</li> </ul>	<ul style="list-style-type: none"> <li>- Clean the adapter or replace if damaged (➤ p. 13-16)</li> </ul>
Recharging procedure properly: LED does not light during charging cycle	<ul style="list-style-type: none"> <li>- Instrument or charger failed</li> </ul>	<ul style="list-style-type: none"> <li>- Send the instrument in for repair or replace the charger (➤ p. 25)</li> </ul>
Recharging procedure properly: LED does light up during charging cycle	<ul style="list-style-type: none"> <li>- Battery failed</li> </ul>	<ul style="list-style-type: none"> <li>- Replace the battery (➤ p. 18-22)</li> </ul>

If recommendations in the Trouble Shooting do not solve current problems, contact your authorized dealer or the manufacturer directly.

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## Return for repair

### **Important!**

Transporting of hazardous materials without a permit is a violation of federal law.

BrandTech Scientific, Inc. will not accept instruments that are not appropriately cleaned and decontaminated.

- Therefore, contact BrandTech Scientific, Inc. and obtain return authorisation **before** sending your instrument for service.

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## Warranty

As provided by law, any and all warranties are null and void if the product has been misused, modified or repaired by unauthorized personnel, if the defects are caused by negligence (instruction manual, maintenance) or by normal wear and tear. Use only original manufacturer's accessory/spare parts.

For conditions and extend of warranty refer to our warranty card enclosed with the product.

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**Best.-Nr. 2 0265 06**

## Ordering Data

### **accu-jet® pipette controller with battery charger,**

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**Wandhalter** Best.-Nr. 265 40

## Spare Parts

**Silicone adapter  
with check valve** Cat. No. 265 08

**Adapter support** Cat. No. 265 12

**Spare membrane filter 0.2 µm, sterile**  
10 pc. in blister pack Cat. No. 2 0265 39

**Spare membrane filter 0.2 µm, non-sterile**  
10 pc. in PE-bag Cat. No. 265 35

**Spare membrane filter 3 µm, non-sterile**  
1 pc. in blister pack Cat. No. 260 52

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This sign certifies that the product meets the requirements of the EEC directive and has been tested according to the specified test methods.